**HAL**  
open science

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Temperature (°C)	Gas (sccm)	Setpoint	Readout	Pressures (mBar)	Recipe parameters	Machine parameters
Control mode: <b>POWER control</b>	N2: 0.0	0.0	0.0	Full Range: <b>NA</b>	Recipe time: 0.0	Idle
Setp. [%]: 0.0	O2: 0.0	0.0	0.0	Capa: 1.33E+1	Step time: 0.0	System DN
Power [%]: 0.0	Ar: 0.0	0.0	0.0		Step number: 0	Cooling: <b>Not cooling</b>
TC1: <b>Absent</b>					Step comment:	Process: <b>Idle</b>
Pyro1: 3.0					Condition:	Chamber: <b>Closed</b>
TC2: <b>Absent</b>					Recipe name: Segton 1000 30s	

Last downloaded recipe		Recipe to download	
Name	Segton 1000 30s	Name	<input type="text"/>
Date of last download	07/09/2015	Creation date	
Hour of last download	15:17:56	Modification date	
User name	Adm	User name	
PID table	Segton-FID-SU-200-picots 2mm	PID table	
TC1 calibration table	Default_TC_Table	TC1 calibration table	
TC2 calibration table	Default_TC_Table	TC2 calibration table	
Pyrometer 1 calibration table	CAL_SU_200mm_2014-09-24-picots 2mm	Pyrometer 1 calibration table	
IR Compensation table	532_ASMAS1_EH2000_S115U-1_AAS-Int4	IR Compensation table	
Post recipe action	Purge	Post recipe action	
Furnace configuration	2000	Furnace configuration	
Recipe extracted from : C:\AnnealSys\Historicals\Segton 1000-500 P 30.his			

Machine	Chamber
Start purge	Open chamber
Purge until ATM	

Recipes	<b>Process</b>	Historicals	Configuration	Manual Mode	Diagnostics	Shut down	Adm/Administrator
---------	----------------	-------------	---------------	-------------	-------------	-----------	-------------------

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Temperature [°C]		Gas (sccm)	Setpoint	Readout	Pressures (mBar)		Recipe parameters		Machine parameters	
Control mode	POWER control	N2	0.0	0.0	Full Range	86	Recipe time	0.0	Idle	
Setp. [%]	0.0	O2	0.0	0.0	Capa	1.33E+1	Step time	0.0	System ON	
Power [%]	0.0	Ar	0.0	0.0			Step number	0	Cooling	Not cooling
TC1	Absent					Step comment			Process	Idle
Pyro1	1.					Condition			Chamber	Closed
TC2	Absent					Recipe name	Segton 1000 30s			

Last downloaded recipe		Recipe to download	
Name	Segton 1000 30s	Name	Segton 1000 30s
Date of last download	07/09/2015	Creation date	07/06/2011 - 14:49:25
Hour of last download	15:17:56	Modification date	07/09/2015 - 15:19:33
User name	Adm	User name	Adm
PID table	Segton-PID-SU-200-picots 2mm	PID table	Segton-PID-SU-200-picots 2mm
TC1 calibration table	Default_TC_Table	TC1 calibration table	Default_TC_Table
TC2 calibration table	Default_TC_Table	TC2 calibration table	Default_TC_Table
Pyrometer 1 calibration table	CAL_SU_200mm_2014-0924-picots 2mm	Pyrometer 1 calibration table	CAL_SU_200mm_2014-0924-picots 2mm
IR Compensation table	9323_ASMAS1EH2000_S115U-1AAS-Int4	IR Compensation table	9323_ASMAS1EH2000_S115U-LAAS-Int4
Post recipe action	Purge	Post recipe action	Purge
Furnace configuration	2000	Furnace configuration	2000

Machine	Chamber
Start purge	Open chamber
Purge until ATM	

Recipes	Process	Historicals	Configuration	Manual Mode	Diagnostics	Shut down	Adm/Administrator Software version: 3.7.0 PCD Version: 3.7.0 09/10/2015 - 09:55:34
---------	---------	-------------	---------------	-------------	-------------	-----------	---

Figure 6: Validation d'une recette

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Temperature [°C]		Gas (sccm)	Setpoint	Readout	Pressures (mBar)		Recipe parameters		Machine parameters	
Control mode	POWER control	N2	0.0	0.0	Full Range	NA	Recipe time	0.0	Idle	
Setp. [%]	0.0	O2	0.0	0.0	Capa	1.33E+1	Step time	0.0	System ON	
Power [%]	0.0	Ar	0.0	0.0			Step number	0	Cooling	Not cooling
TC1	Absent					Step comment			Process	Finished
Pyro1	6.1					Condition			Chamber	Opened
TC2	Absent					Recipe name	Segton 1000+700			

**AnnealSys Control Software**

**Warning**

Manual mode allows direct control of the heating, vacuum and gas equipment. This mode is dedicated to maintenance operation and some safety interlocks are not activated in this mode. Manual mode operations must only be carried out by persons who have been trained by AnnealSys or its representative field service engineers and who have the necessary experience of the system, the processes, the gases and the related risks. Special precautions must be taken in order to avoid dangerous gas mixtures inside the process chamber, or heating with potential explosive gas mixture inside the process

Ok

Software version: 3.7.0  
PCD Version: 3.7.0

Recipes	Process	Historicals	Configuration	Manual Mode	Diagnostics	Shut down	Adm/Administrator Software version: 3.7.0 PCD Version: 3.7.0 11/24/2015 - 15:07:36
---------	---------	-------------	---------------	-------------	-------------	-----------	---

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Recipes	<b>Process</b>	Historicals	Configuration	Manual Mode	Diagnostics	Shut down	Adm/Administrator Software version : 3.7.0 PCD Version : 3.7.0 11/24/2015 - 15:03:51
---------	----------------	-------------	---------------	-------------	-------------	-----------	---

Temperature (°C)	Gas (sccm)	Setpoint	Readout	Pressures (mBar)	Recipe parameters	Machine parameters
Control mode: POWER control	N2	0.0	0.0	Full Range NA	Recipe time: 0.0	Idle
Setp. (%): 0.0	O2	0.0	0.0	Capa: 1.33E+1	Step time: 0.0	System ON
Power (%): 0.0	Ar	0.0	0.0		Step number: 0	Cooling: Not cooling
TC1: Absent					Step comment:	Process: Finished
Pyro1: 1.					Condition:	Chamber: Opened
TC2: Absent					Recipe name: Segton 1000+700	

Last downloaded recipe		Recipe to download	
Name	Segton 1000+700	Name	<input type="text"/>
Date of last download	23/11/2015	Creation date	
Hour of last download	13:42:43	Modification date	
User name	Adm	User name	
PID table	Segton-PID-SU-200-picots 2mm	PID table	
TC1 calibration table	Default_TC_Table	TC1 calibration table	

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Temperature (°C)	Gas (sccm)	Setpoint	Readout	Pressures (mBar)	Recipe parameters	Machine parameters
Control mode: POWER control	N2	0.0	0.0	Full Range NA	Recipe time: 0.0	Idle
Setp. (%): 0.0	O2	0.0	0.0	Capa: 1.33E+1	Step time: 0.0	System ON
Power (%): 0.0	Ar	0.0	0.0		Step number: 0	Cooling: Not cooling
TC1: Absent					Step comment:	Process: Finished
Pyro1: 3.0					Condition:	Chamber: Opened
TC2: Absent					Recipe name: Segton 1000+700	

Last downloaded recipe		Recipe to download	
Name	Segton 1000+700	Name	Segton 1000+700
Date of last download	23/11/2015	Creation date	07/06/2011 - 14:49:25
Hour of last download	13:42:43	Modification date	24/11/2015 - 10:36:24
User name	Adm	User name	Adm
PID table	Segton-PID-SU-200-picots 2mm	PID table	Segton-PID-SU-200-picots 2mm
TC1 calibration table	Default_TC_Table	TC1 calibration table	Default_TC_Table
TC2 calibration table	Default TC Table	TC2 calibration table	Default TC Table

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Temperature [°C]	Gas (sccm)	Setpoint	Readout	Pressures (mBar)	Recipe parameters	Machine parameters
Control mode: POWER control	N2: 0.0	0.0	0.0	Full Range: NA	Recipe time: 0.0	Idle
Setp. [3]: 0.0	O2: 0.0	0.0	0.0	Capa: 1.33E+1	Step time: 0.0	System ON
Power [3]: 0.0	Ar: 0.0	0.0	0.0		Step number: 0	Cooling: Not cooling
TC1: Absent					Step comment:	Process: Finished
Pyro1: 4.6					Condition:	Chamber: Opened
TC2: Absent					Recipe name: Segton 1000+700	

**Recipe header information**

System: AS-MASTER  
 Creation date: 07/06/2011 - 14:49:25  
 Modification date: 24/11/2015 - 10:36:24  
 User Name: Adm

Recipe name: **Segton 1000+700**

TC1 calibration table: Default\_TC\_Table  
 TC 2 calibration table: Default\_TC\_Table  
 Pyrometer 1 calibration table: CAL\_SU\_200mm\_2014-0924-picots  
 PID control parameter table: Segton-PID-SU-200-picots 2mm  
 IR compensation table: 5323\_ASMASTER2000\_S1150-LAAS  
 Furnace configuration: 2000 HT  
 Post recipe action: Purge Clean None

Buttons: New, Browse, Delete, Print, Save as, Save

Bottom navigation: Recipes, Process, Historicals, Configuration, Manual Mode, Diagnostics, Shut down

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**Validation**

Confirmation dialog: Overwrite file C:\AnnealSys\Recipes\Segton 1000+700 already exists. Overwrite it? [Ok] [No]

**Confirmation**

Information dialog: The file has been successfully saved. [Ok]

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